

PAGEr™ Gel Spacer for the XCell SureLock® Chamber

Instructions

PAGEr™ (10 cm x 10 cm) Gels perform effectively in the Novex® XCell SureLock® Mini-Cell when used with the Lonza PAGEr™ Gel Spacer.

To run one gel: Put the gel in the front of the chamber. Use Novex® buffer dam on the back. Place the Lonza spacer between the gel and the buffer core, on the side of the chamber with the gel tension wedge. Lock the gel tension wedge.

To run two gels: Put a gel on each side of the buffer core. Place the Lonza spacer between the gel and the buffer core, on the side of the chamber with the gel tension wedge. Lock the gel tension wedge.

Ordering Information

Catalog No.	Description
59900	PAGEr™ Gel Spacer Spacers sent free of charge.

Related Products

PAGEr™ Gold Precast Gels 10 cm x 10 cm

PAGEr™ EX Precast Gels 10 x 10 cm

PAGEr™ Minigel Chamber Catalog No. 59905.

(see www.lonza.com for complete listing)

Manufactured for Lonza.

XCell SureLock and Novex are trademarks of Life Technologies Corp. All other trademarks herein are marks of the Lonza Group or its affiliates.

© 2012 Lonza Rockland, Inc.
All rights reserved.