

www.lonza.com
U.S. Scientific Support: 800-521-0390
scientific.support@lonza.com
EU/ROW Scientific Support: +49-221-99199-400
scientific.support.eu@lonza.com
Document # INST-12-725 11/11
© 2011 Lonza Walkersville, Inc.

## BioWhittaker<sup>®</sup> UltraCULTURE™ Serum-free Medium

#### Introduction

UltraCULTURE<sup>™</sup> Medium (cat. no. 12-725F) is a serum-free medium designed for the cultivation of a wide variety of mammalian cell types. The medium consists of a DMEM:F-12 base which is supplemented with recombinant human insulin, bovine transferrin and a purified mixture of bovine serum proteins including albumin. The total protein concentration of UltraCULTURE<sup>™</sup> Medium is approximately 3 mg/ml. UltraCULTURE<sup>™</sup> Medium does not contain L-glutamine.

UltraCULTURE™ Medium has been used to grow a variety of cells of primary origin and established cell lines. The medium may be used to support the fusion of cells during the formation of hybridomas. UltraCULTURE™ Medium has been used to grow cells of lymphoid origin including monocyte and macrophage cell lines. Epithelial and fibroblastic cells have been cultivated in UltraCULTURE™ Medium and endothelial cells have been cultivated under reduced serum concentration (5% fetal bovine serum using UltraCULTURE™ Medium as a base formulation).

UltraCULTURE™ Medium has also been used as a complete medium to support the generation of viral particles for use in vaccine production.

UltraCULTURE™ Medium may be supplemented with cryoprotective medium (cat. no. 12-132) to cryopreserve cells in a serum-free environment.

#### Instructions for Use

UltraCULTURE™ Medium is a complete all-purpose serum-free medium which supports the growth of a wide variety of both non-adherent and adherent cell lines. The product is ready-to-use with the simple addition of 5.0 ml of L-glutamine solution (cat. no. 17-605) per 500 ml.

UltraCULTURE  $^{\text{TM}}$  Medium should be stored at  $2^{\circ}$  -  $8^{\circ}$ . For cells in serum-containing medium weaning is NOT required.

Turbidity may develop UltraCULTURE™ Medium. Experiments have determined that the turbidity will not alter the performance of the product.

# Partial List of Cells Cultivated with UltraCULTURE™ Medium:

Ultracolitore Medium.				
Cell line	Source	Cell type		
HEL, N-10	Human	Fetal lung diploid		
		fibroblast		
HeLa	Human	Uterine cancer		
HuL-1,2	Human	Liver (normal)		
HuK-1	Human	Kidney (normal)		
HuS-1AT	Human	Skin		
HEC	Human	Embryonic cancer		
HL-60	Human	Acute promyelocytic leukemia		
Raji	Human	Burkitt's lymphoma		
EB-3	Human	Burkitt's lymphoma		
K-562	Human	Chronic myelocytic leukemia		
HNK	Human	Neonatal kidney (primary)		
HTC29	Human	Colon Cancer		
TT	Human	Medulary thyroid tumor		
MB231	Human	Breast carcinoma		
U138	Human	Glioma		
FM3A	Mouse	Breast cancer		
NS-1	Mouse	Myeloma		
L	Mouse	Subcutaneous		
P388D1	Mouse	Macrophage-like		
P815	Mouse	Mast cell tumor		
T3	Mouse	Pituitary		
B82	Mouse	L cells –connective tissue		
RPL-1	Rat	Peritoneum		
RSP-2	Rat	Spleen		
RLG-1	Rat	Lung		
Lym-1	Rat	Lymph node		
RCR-1	Rat	Brain		
235-1, MNQ	Rat	Pituitary		
GC, GH3	Rat	Pituitary		
CA77	Rat	Medulary thyroid tumor		
Rat-1	Rat	Fibroblast		
JTC-12	Monkey	Kidney		
COS1, COS7	African green monkey	SV40 transformed kidney		



## Storage

2℃ to 8℃

#### **Product Use Statement**

### THESE PRODUCTS ARE FOR RESEARCH USE

**ONLY.** Not approved for human or veterinary use, for application to humans or animals, or for use in clinical or *in vitro* procedures.

## **Ordering Information**

Catalog	Description	Size
number		
12-725F	UltraCULTURE <sup>™</sup> Medium	500 ml
	without L-glutamine	