


Amaxa[®] Human Dermal Fibroblasts Nucleofector[®] Kit


For Normal Human Dermal Fibroblasts – Neonatal (NHDF-Neo)

Validated to work with Clonetics[®] NHDF-Neo [Lonza; Cat. No. CC-2509]; adherent fibroblastoid cells

Example for Nucleofection[®] of NHDF-Neo

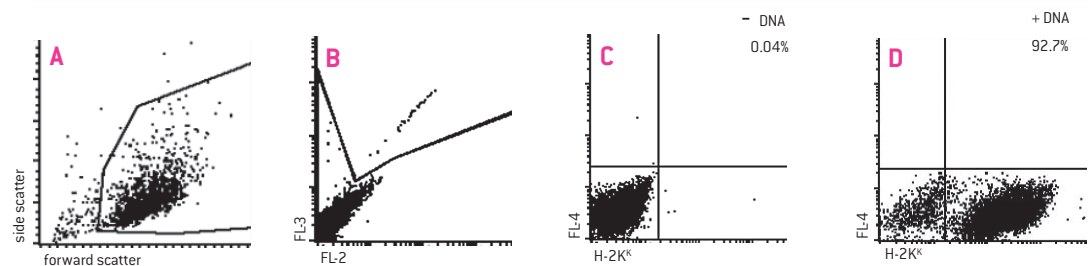

NHDF-Neo were transfected using the Human Dermal Fibroblast Nucleofector[®] Kit, program U-020 and a plasmid encoding the mouse MHC class I heavy chain molecule, H-2K^k. 5 hours post Nucleofection[®] the cells were stained with a Cy5-coupled antibody directed against H-2K^k and were analyzed by flow cytometry. NHDF-Neo were gated according to forward/side scatter (A). Dead cells were excluded by staining with propidium iodide and gating (B). H-2K^k expression of NHDF-Neo is shown after Nucleofection[®] without (C) and with plasmid DNA (D).

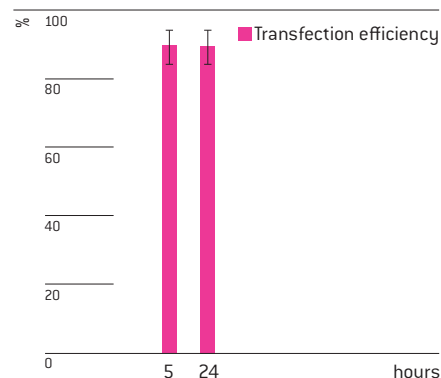

Transfection efficiencies of NHDF-Neonatal 5 and 24 hours post Nucleofection[®]. Cells were transfected using program U-020 and a plasmid encoding the mouse MHC class I heavy chain molecule H-2K^k. Viability is usually around 85 – 90%.

Product Description

Cat. No.	VPD-1001
Size (Reactions)	25
Human Dermal Fibroblast Nucleofector [®] Solution	2.25 ml (2.05 ml + 10% overfill)
Supplement	0.5 ml (0.45 ml + 10% overfill)
pmaxGFP [®] Vector (0.5 µg/µl in 10 mM Tris pH 8.0)	30 µg
Certified cuvettes	25
Plastic pipettes	25
Storage and stability	Store Nucleofector [®] Solution, Supplement and pmaxGFP [®] Vector at 4°C. For long-term storage, pmaxGFP [®] Vector is ideally stored at -20°C. The expiration date is printed on the solution box. Once the Nucleofector [®] Supplement is added to the Nucleofector [®] Solution it is stable for three months at 4°C.

Required Material

Note Please make sure that the entire supplement is added to the Nucleofector® Solution. The ratio of Nucleofector® Solution to supplement is 4.5 : 1. For a single reaction use 82 µl of Nucleofector® Solution plus 18 µl of supplement to make 100 µl of total reaction volume.

- Nucleofector® Device
- Supplemented Nucleofector® Solution at room temperature
- Supplied certified cuvettes
- Supplied plastic pipettes
- Supplied pmaxGFP® Vector
- Substrate of interest, highly purified, preferably by using endotoxin free Kits; A260 : A280 ratio should be at least 1.8
- 6-well culture dish
- **For trypsinization:** Reagent Pack™ Subculture Reagent Kit containing Trypsin/EDTA, HEPES Buffered Saline Solution (HBSS) and Trypsin Neutralizing Solution (TNS) [Lonza, Cat. No. CC-5034]. Alternatively if cells hardly detach: Trypsin 0.5 % – EDTA 0.2%
- **Culture medium:** FGM®-2 BulletKit [Lonza; Cat. No. CC-3132]. **We recommend storing 40 ml aliquots of the prepared medium at -80°C. Do not use medium stored at 4°C for more than two or three days, as this may lead to reduced cell viability**
- Prewarm appropriate volume of culture medium to 37°C (1.5 ml per sample)
- Appropriate number of cells (0.5 – 1 x 10⁶ cells per sample)
Minimal cell number: 2 x 10⁵ cells (a lower cell number may lead to a major increase in cell mortality)
Maximum cell number: 2 x 10⁶

1. Pre Nucleofection®

Note Transfection results may be donor-dependent.

Cell culture recommendations

- 1.1 Seeding conditions: at least 5 x 10⁴ cells/25cm² flask
- 1.2 Replace media 2 – 3 times per week; 2 – 3 ml media per 25 cm² flask
- 1.3 Cells should be passaged after reaching 90% confluency
- 1.4 Cells should be preferably passaged 2 days before Nucleofection®
- 1.5 Do not use cells after passage number 15 as this may result in substantially lower gene transfer efficiency and viability
- 1.6 Optimal confluency before Nucleofection® 90%

Trypsinization

- 1.7 Remove media from the cultured cells and wash cells once with HBSS; use at least same volume of HBSS as culture media
- 1.8 For harvesting, incubate the cells 1 – 3 minutes at 37°C with recommended volume of indicated trypsinization reagent (please see required material)
- 1.9 Neutralize trypsinization reaction with TNS once the majority of the cells (>90%) have been detached

2. Nucleofection®

One Nucleofection® Sample contains

0.5 – 1 x 10 ⁶ cells
1 – 5 µg plasmid DNA (in 1 – 5 µl H ₂ O or TE) or 2 µg pmaxGFP® Vector or 30 – 300 nM siRNA (3 – 30 pmol/sample)
100 µl Human Dermal Fibroblast Nucleofector® Solution

- 2.1 Please make sure that the entire supplement is added to the Nucleofector® Solution
- 2.2 Prepare 6-well plates by filling appropriate number of wells with 1.5 ml of supplemented culture media and pre-incubate/equilibrate plates in a humidified 37°C/5% CO₂ incubator
- 2.3 Harvest the cells by trypsinization (please see 1.7 – 1.9)
- 2.4 Count an aliquot of the trypsinized cells and determine cell density
- 2.5 Centrifuge the required number of cells (0.5 – 1 x 10⁶ cells per sample) at 200xg for 10 minutes at room temperature
- 2.6 Resuspend the cell pellet carefully in 100 µl room temperature Nucleofector® Solution per sample

Note Avoid storing the cell suspension for longer than 15 minutes in Nucleofector® Solution as this reduces cell viability and transfection efficiency.

- 2.7 Combine 100 µl of cell suspension with 1 – 5 µg DNA, 2 µg pmaxGFP® Vector or 30 nM – 300 nM siRNA (3 – 30 pmol/sample) or other substrates
- 2.8 Transfer cell/DNA suspension into certified cuvette (sample must cover the bottom of the cuvette without air bubbles. Close the cuvette with the cap)
- 2.9 Select the appropriate Nucleofector® Program U-020 (U-20 for Nucleofector® I Device)
- 2.10 Insert the cuvette with cell/DNA suspension into the Nucleofector® Cuvette Holder and apply the selected program
- 2.11 Take the cuvette out of the holder once the program is finished
- 2.12 Add ~500 µl of the pre-equilibrated culture media to the cuvette and gently transfer the sample immediately into the 6-well plate (final volume 1.5 ml media per well/sample). Use the supplied pipettes and avoid repeated aspiration of the sample

3. Post Nucleofection®

- 3.1 Incubate the cells in a humidified 37°C/5% CO₂ incubator until analysis. Gene expression is often detectable already after 4 – 5 hours

Additional Information

For an up-to-date list of all Nucleofector® References, please refer to:
www.lonza.com/nucleofection-citations

For more technical assistance, contact our Scientific Support Team:

USA/Canada

Phone: 800 521 0390 (toll-free)

Fax: 301 845 8338

E-mail: scientific.support@lonza.com

Europe and Rest of World

Phone: +49 221 99199 400

Fax: +49 221 99199 499

E-mail: scientific.support.eu@lonza.com

References

1. Frolov A, Zielinski et al, J Biol Chem. 2003;278(28):25517-25.
2. Javelaud D et al, J Biol Chem. 2003;278(27):24624-8.
3. Neff L et al, J Biol Chem. 2003;278(30):27721-8.
4. Ohtani N et al, J Cell Biol. 2003;162(2):173-83.

Lonza Cologne AG
50829 Cologne, Germany

Please note that the Amaxa® Nucleofector® Technology is not intended to be used for diagnostic purposes or for testing or treatment in humans.

The Nucleofector® Technology, comprising Nucleofection® Process, Nucleofector® Device, Nucleofector® Solutions, Nucleofector® 96-well Shuttle® System and 96-well Nucleocuvette® plates and modules is covered by patent and/or patent-pending rights owned by Lonza Cologne AG.

Amaxa, Nucleofector, Nucleofection, maxGFP, Clonetics, Reagent Pack, FGM and BulletKit are either registered trademarks or trademarks of the Lonza Group or its affiliates.

Other product and company names mentioned herein are the trademarks of their respective owners.

This kit contains a proprietary nucleic acid coding for a proprietary copepod fluorescent protein intended to be used as a positive control with this Lonza product only. Any use of the proprietary nucleic acid or protein other than as a positive control with this Lonza product is strictly prohibited. USE IN ANY OTHER APPLICATION REQUIRES A LICENSE FROM EVROGEN. To obtain such a license, please contact Evrogen at license@evrogen.com.

The CMV promoter is covered under U.S. Patents 5,168,062 and 5,385,839 and its use is permitted for research purposes only. Any other use of the CMV promoter requires a license from the University of Iowa Research Foundation, 214 Technology Innovation Center, Iowa City, IA 52242.

The use of this product in conjunction with materials or methods of third parties may require a license by a third party. User shall be fully responsible for determining whether and from which third party it requires such license and for the obtaining of such license.

No statement is intended or should be construed as a recommendation to infringe any existing patent.

© Copyright 2009, Lonza Cologne AG. All rights reserved DPD-1001 09/09